


FISH PROBES - OACP BUFFERS COMPATIBILITY

READY TO USE FISH
PROBES

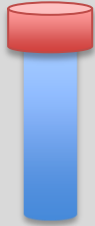

The Smart-ISH and Rapid-ISH hybridization buffers are compatible with both ready to use probes and those that need to be reconstituted. Just add 5ul of the OaCP Buffer to the probe prepared as usual.
www.oacp.it- ask@oacp.it



PROBES THAT NEED
TO BE
PREPARED/DILUITED



RECONSTITUTE
AS FROM PROBE
INSTRUCTIONS *



ADD 1-5UL PROBE
ON
THE SLIDE



ADD 5UL OF
OACP BUFFER
ON THE
SLIDE, MIX



DENATURATION
AND
HYBRIDATION
STEPS

*The OaCP Buffers **do not replace** the elution buffer already present in the package.